



Removing the discharge protector

Prior to first usage please proceed as follows:

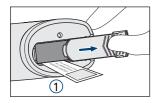

Slide the battery compartment lid open.

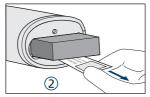


Remove the battery and discharge protection.




Install the battery and close the battery compartment lid.

Battery status display


> Ready for operation

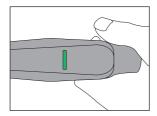

Solid green light: The rechargeable battery is fully charged.

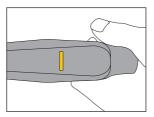


Solid yellow light: The rechargeable battery is half-charged.

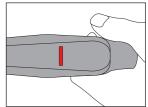

Solid red light: The rechargeable battery is almost empty.

> Empty battery

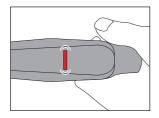

Flashing red light: The rechargeable battery needs to be charged.


Loading battery






Battery status display


> Charge battery

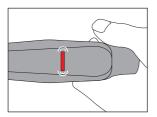

Flashing red light: The rechargeable battery is <5% charged.

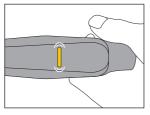

Flashing yellow light: The rechargeable battery is 5-60% charged.

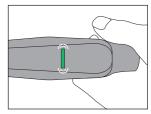


Flashing green light: The rechargeable battery is 60-99% charged.

Speed control


The liquid aspiration speed is regulated continuously by controlling the pressure on the control buttons, i.e. how far in the button is pressed.

> Slow aspiration

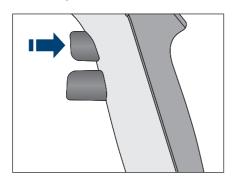

To slowly aspirate or dispense liquid, press the corresponding control button lightly.

> Quick aspiration

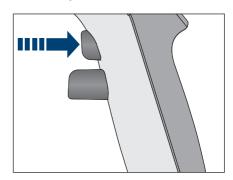

To quickly aspirate or dispense liquid, press the corresponding control button firmly.