



## Autoclaving Instructions for TipOne<sup>®</sup> Products

1. Place racked tips in the autoclave. If desired, the hinged lid may be propped open using the tab on the front of the lid.
2. Autoclave at 121 °C (252 °F) under a pressure of 15 psi for 15 minutes.
3. Allow the tips to dry thoroughly before use.

**Note:**

Remove the shrink-wrap from stacked tips before autoclaving.

Tips in reload packages must be placed inside the hinged racks before autoclaving.

Excessive heat and/or prolonged time in the autoclave may cause warping of both the tips and the rack.