



LASER CRYO-BABIES[®] AND CRYOTAGS[®]

Labels for Cryogenic Storage*


DURABLE - Chemically inert, polyolefin labels resist most organic solvents and caustic agents. They adhere to most plastics†, glass, and metals without cracking, peeling, or degrading.

TEMPERATURE-RESISTANT - From -196°C to 100°C, Cryo-Babies and Cryo-Tags withstand boiling water baths, moderate oven temperatures (150°C), autoclaving and extreme freezer temperatures (liquid and vapor phase nitrogen).

SPECIALLY DESIGNED - Label sheets are specially top-coated to readily accept laser printing. Heat-resistant, stay-flat sheets will not jam in laser printers.

IDEALLY SIZED - Will fit a wide variety of cryovials, microtubes, storage boxes and other containers.

COLOR CODING - Each label size is available in white, blue, green, orange, red and yellow.

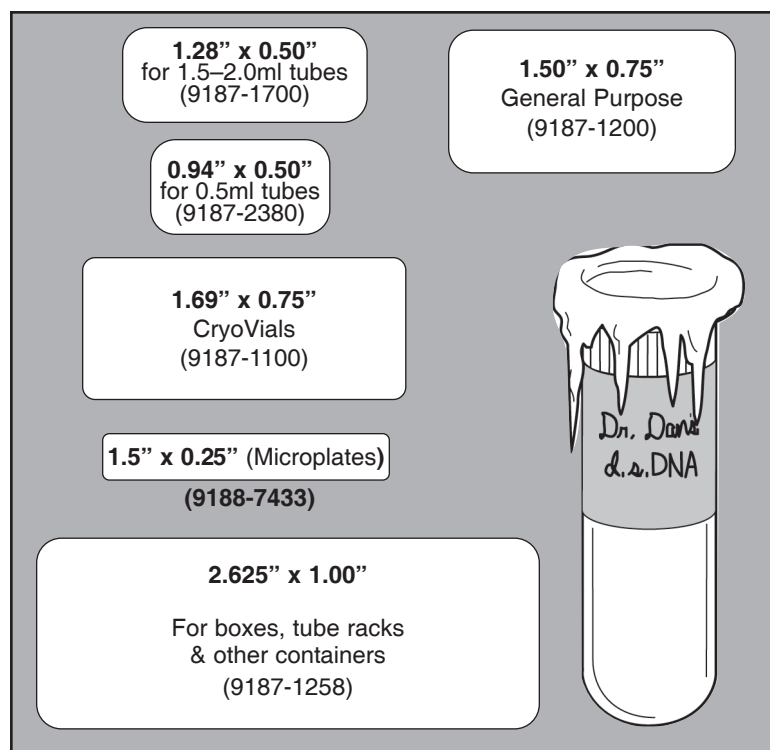

CAT. NO.	COLOR	LABEL SIZE	LABELS PER SHT/PK	PRICE
9187-1100	Specify	1.69" x 0.75" / 43mm x 19mm	52 / 1,040	60.00
9187-1200	Specify	1.50" x 0.75" / 38mm x 19mm	60 / 1,200	60.00
9187-1700	Specify	1.28" x 0.50" / 33mm x 13mm	85 / 1,700	60.00
9187-2380	Specify	0.94" x 0.50" / 24mm x 13mm	119 / 2,380	60.00
9188-7433	White	1.50" x 0.25" / 38mm x 6mm	156 / 3,120	60.00
9187-1258	White	2.625" x 1.00" / 67mm x 25mm	30 / 600	60.00
9187-1108	Assorted*	1.69" x 0.75" / 43mm x 19mm	52 / 1,040	60.00
9187-1208	Assorted*	1.50" x 0.75" / 38mm x 19mm	60 / 1,200	60.00
9187-1708	Assorted*	1.28" x 0.50" / 33mm x 13mm	85 / 1,700	60.00
9187-2388	Assorted*	0.94" x 0.50" / 24mm x 13mm	119 / 2,380	60.00
9188-7408	Assorted*	1.50" x 0.25" / 38mm x 6mm	156 / 3,120	60.00
9187-2508	Assorted*	2.625" x 1.00" / 67mm x 25mm	30 / 600	60.00

For a Color (*other than white*) replace the last digit of the catalog number with: 1=Blue, 2=Green, 3=Orange, 4=Red, 6=Yellow (Example: 9187-1701 for Blue)

*Assorted Packs include 4 sheets of each color: Blue, Green, Orange, Red and Yellow

ORDER INFORMATION

Online: www.usascientific.com

Phone: 1-800-522-8477
1-352-237-6288

Fax: 1-352-351-2057

Mail: USA Scientific, Inc.
PO Box 3565
Ocala, FL 34478-3565

† NOTE: Siliconized surfaces should be avoided. Apply to clean, dry tubes at room temperature. Tubes made of 100% virgin polypropylene are recommended.

*patented

See reverse side for formatting instructions in Microsoft Word[®] →



PO Box 3565 Ocala, FL 34478-3565
 Phone:1-800-522-8477 or 1-352-237-6288
 Fax:1-352-351-2057 Web: www.usascientific.com

FORMATTING INSTRUCTIONS FOR
LASER CRYO-BABIES®
 and **CRYO-TAGS®**
 IN MICROSOFT WORD®

1. Open Word
2. Click on Tools > Letters & Mailings (if you have Office XP. Office 2000 users skip to step 4.)
3. Click on Envelopes and Labels > Labels tab
4. Click on Options
5. Under Printer Information select the Laser option
6. Under Tray, select Manual Feed
7. Click on New Label to customize your label
8. Type in a name for the label (e.g. Catalog Number)
9. Insert the Label Details which can be found at the bottom of this page and click OK
10. Select the label you made in the Label Options window and click OK
11. In the Envelopes & Labels window, type your information in the area provided and click on New Document
12. When the document appears on the screen, click Edit > Select All
13. When the whole document is highlighted, choose an appropriate font and font size* and center the type using the Center command located in the toolbar

If the text in either the rows or columns do not align, continue with Steps 14 -16 otherwise skip to Step 17.

- | | |
|-----|--|
| 14. | After centering the text, display the Tables and Borders Toolbar by clicking View>Toolbars>Tables and Borders |
| 15. | Highlight the top half of the sheet and click on the Align Center icon in the Tables and Borders Toolbar |
| 16. | Next, highlight the bottom half of the sheet and click the Align Top Center, Left or Right (depending on where the text needs to be aligned) icon in the Tables and Borders Toolbar. |
17. To determine if your text will fall within the boundaries of the label, have the Gridlines showing
 18. To make the gridlines appear on the screen, click on Table > Show Gridlines
 19. Before printing a sheet of labels, print a test page on blank paper and line it up with a sheet of labels to ensure the text lines up correctly

Notes:

- Pass the laser sheet through the printer only once
 - Manually feed the laser sheets one at a time
 - Print your labels on a blank sheet of paper and check the positioning on a blank sheet of labels
 - Minor adjustments may be required depending on your printer and/or software you are using
- * Suggested Font Sizes: 8 pt. for smaller labels or 10-12 pt. for larger labels

Catalog No.	9187-1700	9187-1200	9187-2380	9187-1100	9187-1258	9188-7433
Label Size	1.28 x 0.50"	1.50 x 0.75"	0.94 x 0.50"	1.69 x 0.75"	2.625 x 1.00"	1.50 x 0.25"
Top Margin	0.24"	0.31"	0.24"	0.63"	0.50"	0.62"
Side Margin	0.77"	0.25"	0.56"	0.70"	0.19"	0.50"
Vert. Pitch	0.63"	0.88"	0.63"	0.75"	1.00"	0.25"
Horiz. Pitch	1.40"	1.63"	1.07"	1.80"	2.75"	2.00"
Label Height	0.50"	0.75"	0.50"	0.75"	1.00"	0.25"
Label Width	1.28"	1.50"	0.94"	1.69"	2.625"	1.50"
No. Across	5	5	7	4	3	4
No. Down	17	12	17	13	10	39
Page Size	8.5" x 11"	8.5" x 11"	8.5" x 11"	8.5" x 11"	8.5" x 11"	8.5" x 11"

See reverse side for more information on Cryo-Babies® and Cryo-Tags® →