


## Quick Adjustment Guide for the USA Scientific Repeating Pipet

Volume Selection Dial					Tip Choice
1	2	3	4	5	
Sample Volume					
10 $\mu$ l	20 $\mu$ l	30 $\mu$ l	40 $\mu$ l	50 $\mu$ l	0.5 ml
25 $\mu$ l	50 $\mu$ l	75 $\mu$ l	100 $\mu$ l	125 $\mu$ l	1.25 ml
50 $\mu$ l	100 $\mu$ l	150 $\mu$ l	200 $\mu$ l	250 $\mu$ l	2.5 ml
100 $\mu$ l	200 $\mu$ l	300 $\mu$ l	400 $\mu$ l	500 $\mu$ l	5.0 ml
250 $\mu$ l	500 $\mu$ l	750 $\mu$ l	1.0 ml	1.25 ml	12.5 ml
500 $\mu$ l	1.0 ml	1.5 ml	2.0 ml	2.5 ml	25 ml
1.0 ml	2.0 ml	3.0 ml	4.0 ml	5.0 ml	50 ml
Number of Dispensing Steps					 <a href="http://www.usascientific.com">www.usascientific.com</a>
48	23	15	11	8	